

How to Use the Victor Nivo Plate Reader

1. Turn on the machine and turn on the computer
2. Login to the Victor Nivo computer
3. Open Google Chrome
4. Type in the following code into the navigation bar: **192.168.0.2**
5. Select **Ok** (no need for a username or password)
6. Look for quick guide next to computer in order to properly set up your protocol
7. To analyze data: The software is called MyAssays and is located on the desktop
8. When you are done using the plate reader, **Shut Down** computer and **Turn Off** plate reader (do not just log out—this helps keep our Victor Nivo happy 😊)

If you have any questions please ask a technician in person or email us at ric_tech@byu.edu